

(A) hPSCs cultured in TeSR™-AOF demonstrate a higher plating efficiency compared to hPSCs maintained in low-protein medium (TeSR™-E8™). Plating efficiency is calculated by seeding a known number of aggregates and comparing to the number of established colonies on day 7. (B) hPSCs maintained in TeSR™-AOF exhibit a higher average fold expansion per passage compared to TeSR™-E8™. (C) hPSCs cultured in TeSR™-AOF demonstrate consistent expansion and minimal variability across all cell lines assessed. Cumulative fold expansion was measured from passage 1 to 5. Data represented as mean plating efficiency or fold expansion across 10 passages \pm SD. MG = Matrigel®; VN = Vitronectin XF™.

FEATURE 3 Stabilized Components

Stabilized components, including FGF2, support high cell quality while allowing for restricted feeding schedules. This means that, compared to following a daily feeding schedule, you can save time and media while still obtaining large numbers of high-quality cells.

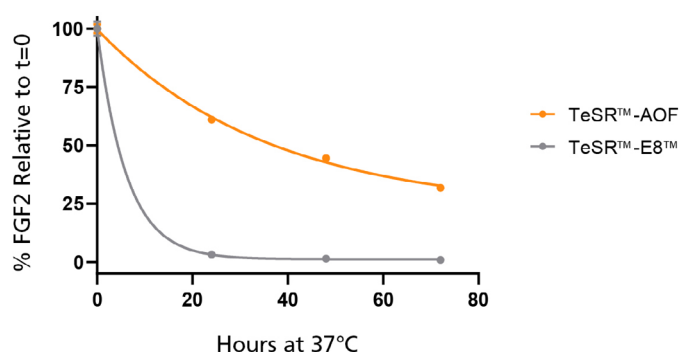


Figure 3. Native FGF Levels Are Stabilized at 37°C in TeSR™-AOF, Allowing Restricted Feeding Schedules

TeSR™-AOF and TeSR™-E8™ were incubated at 37°C for 24, 48, and 72 hours. FGF2 levels were measured by Meso Scale Discovery (MSD) immunoassay; data was normalized to $t = 0$ levels for TeSR™-E8™ and TeSR™-AOF, respectively. FGF2 levels in TeSR™-AOF decline much more slowly than in TeSR™-E8™, with $36.7 \pm 5.61\%$ of $t = 0$ levels at 72 hours when incubated at 37°C. Data representative of $n = 3$ biological replicates \pm SD.

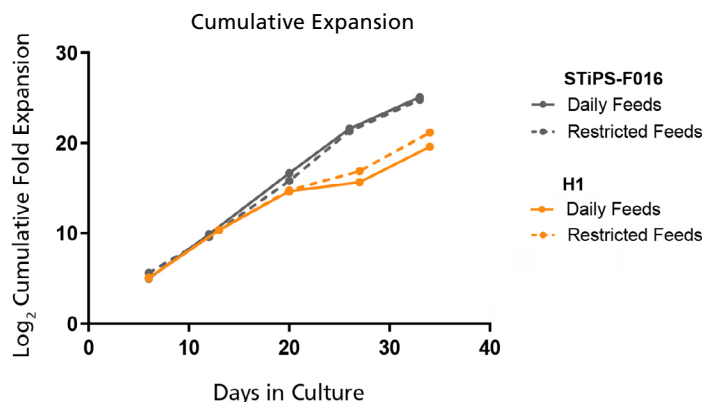


Figure 4. hPSCs Maintained in TeSR™-AOF with Daily and Restricted Feed Schedules Have Comparable Expansion Rates

hPSCs were maintained on Vitronectin XF™ for five passages. At the end of each passage, cell counts were obtained using the Nucleocounter® NC-200 ChemoMetec automated cell counter to count DAPI-stained nuclei. The \log_2 transformed cumulative fold expansion was plotted against time in culture (days).

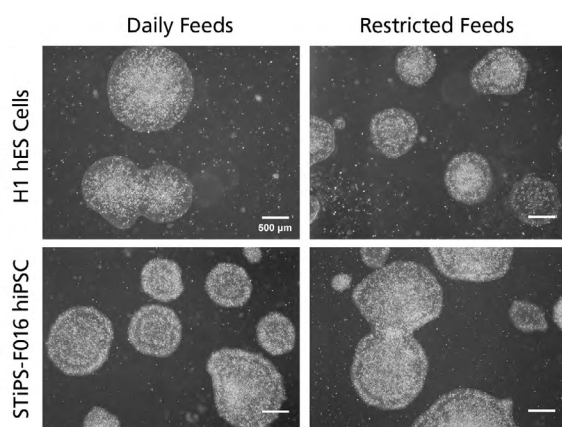


Figure 5. hPSCs Maintained in TeSR™-AOF with Daily and Restricted Feed Schedules Exhibit Comparable Colony Morphology

hPSCs were maintained on Vitronectin XF™ for five passages. Phase-contrast images were taken on day 7 after seeding. For restricted feeds (right panel), hPSCs were fed with a double volume (4 mL) of medium on day 2 after passage, followed by two consecutive skipped days of feeds, with a final single-volume feed (2 mL) on day 5, prior to passaging on day 6 or 7.

FEATURE 4 A Smoother Path to the Clinic

To support cell and gene therapy research, TeSR™-AOF is manufactured under relevant cGMPs with enhanced documentation, including an FDA Master File. To learn more about how cGMP-manufactured TeSR™-AOF can support your clinical applications, visit www.stemcell.com/why-tesr-aof. There, you'll find scientific posters and talks, answers to some frequently asked questions, and a form where you can sign up to request a sample.

For more information on how we can support your regulatory needs, including navigating requirements for using TeSR™-AOF in your cell therapy applications, visit www.stemcell.com/regulatory-support or contact your local STEMCELL representative.

Copyright © 2022 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, are trademarks of STEMCELL Technologies Canada Inc. TeSR, E8, and mTeSR are trademarks of WARF. Vitronectin XF is a trademark of, and developed and manufactured by Nucleus Biologics. Matrigel is a registered trademark of Corning, Inc. Nucleocounter is a trademark of ChemoMetec A/S. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.

PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED. FOR ADDITIONAL INFORMATION ON QUALITY AT STEMCELL, REFER TO WWW.STEMCELL.COM/COMPLIANCE.