

ReproTeSR™

Reprogramming medium for human iPS cell induction

Catalog #05926

500 mL Kit



Scientists Helping Scientists™ | WWW.STEMCELL.COM

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

Product Description

ReproTeSR™ is a complete and xeno-free reprogramming medium developed for generating human induced pluripotent stem (iPS) cells from fibroblasts, urine-derived cells, and blood-derived cells such as CD34+ or erythroid precursor cells under feeder-free conditions.

This medium is recommended for use during the induction phase of reprogramming and may be used with Corning® Matrigel® hESC-Qualified Matrix (Corning Catalog #354277).

Easy to identify and select colonies

- Pre-screened components ensure high quality of iPS cell colony morphology for improved manual selection

Rapid subcloning

- Large colony size and reduced differentiation enable rapid establishment of homogeneous iPS cell cultures

Reproducible efficiency

- Feeder-free, defined formulation facilitates reproducibly efficient human iPS cell generation

Product Information

The following components are sold as a complete kit (Catalog #05926) and are not available for individual sale.

COMPONENT NAME	COMPONENT #	SIZE	STORAGE	SHELF LIFE
TeSR™-E7™/ReproTeSR™ Basal Medium	05919	480 mL	Store at 2 - 8°C.	Stable until expiry date (EXP) on label.
ReproTeSR™ 25X Supplement*	05928	20 mL	Store at -20°C.	Stable until expiry date (EXP) on label.

*This product contains material derived from human plasma. Donors have been tested and found negative for HIV-1 and -2, hepatitis B, and hepatitis C prior to donation. However, this product should be considered potentially infectious and treated in accordance with universal handling precautions.

Preparation of Complete ReproTeSR™ Medium

Use sterile technique when preparing complete ReproTeSR™ medium (Basal Medium + 25X Supplement).

The following example is for preparing 500 mL of ReproTeSR™ medium. If preparing other volumes, adjust accordingly.

1. Thaw 25X Supplement at room temperature (15 - 25°C) or at 2 - 8°C.

NOTE: Once thawed, use immediately or aliquot and store at -20°C. Do not exceed the expiry date as indicated on the label. After thawing the aliquots, use immediately. Do not re-freeze.

2. Add 20 mL of 25X Supplement to 480 mL of Basal Medium. Mix thoroughly.

NOTE: If prepared using sterile technique, complete ReproTeSR™ medium is ready for use and does not require filtering. If not used immediately, store complete ReproTeSR™ medium at 2 - 8°C for up to 2 weeks. Alternatively, aliquot and store at -20°C for up to 1 month. Do not exceed the shelf life of the individual components. After thawing complete medium, use immediately or store at 2 - 8°C for up to 2 weeks. Do not re-freeze.

Directions for Use

Protocols for reprogramming various cell types are described in the documents listed below, available at www.stemcell.com or contact us to request a copy.

Erythroid progenitor cells

Technical Manual for Erythroid Progenitor Reprogramming Kit (Catalog #05924)

CD34+ progenitor cells

Technical Manual for CD34+ Progenitor Reprogramming Kit (Catalog #05925)

Fibroblasts

Product Information Sheet for ReproRNA™-OKSGM Kit (Catalog #05930)



This product was developed under license to intellectual property owned by WiCell™ Research Institute.

This product is sold for research use only (whether the buyer is an academic or for-profit entity) under a non-transferable, limited-use license. Purchase of this product does not include the right to sell, use or otherwise transfer this product for commercial purposes (i.e., any activity undertaken for consideration, such as use of this product for manufacturing, or resale of this product or any materials made using this product, or use of this product or any materials made using this product to provide services) or clinical use (i.e., administration of this product or any material using this product to humans) or the right to implant any material made using this product into an animal by, or in collaboration with, a for-profit entity, for purposes other than basic pre-clinical research applications (including without limitation teratoma assays) to validate the function of the cells. Purchasers who do not agree to the terms and conditions set forth above should return the product in acceptable conditions to the seller for a refund.

PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED. FOR ADDITIONAL INFORMATION ON QUALITY AT STEMCELL, REFER TO WWW.STEMCELL.COM/COMPLIANCE.

Copyright © 2021 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, ReproRNA, and ReproTeSR are trademarks of STEMCELL Technologies Canada Inc. TeSR, E7, and E8 are trademarks of WARF. Corning and Matrigel are registered trademarks of Corning Incorporated. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.