

Easy 250 EasySep™ Magnet

Magnet for column-free immunomagnetic separation

Catalog #100-0821

Catalog #700-0320

Document #10000012623 | Version 00



Scientists Helping Scientists™ | www.stemcell.com

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

Product Description

The Easy 250 EasySep™ Magnet is an anodized aluminum body housing rare earth magnetic elements.

The Easy 250 EasySep™ Magnet is designed for cell separation procedures using EasySep™ reagents. The Easy 250 EasySep™ Magnet generates a high-gradient magnetic field in the interior cavity that is strong enough to separate cells labeled with magnetic EasySep™ particles, without the use of columns. This magnet is designed to hold a standard T-75 cm² flask (i.e. Corning Catalog #353135).

Properties

Storage: Store at 15 - 25°C.

Shelf Life: Not applicable

STRONG MAGNETIC FIELD: Keep away from pacemakers, magnets, computer disks, watches, and other objects that respond to magnetic fields.

Handling / Directions For Use

The Easy 250 EasySep™ Magnet is only for use with certain EasySep™ kits. For more information, refer to the appropriate EasySep™ Product Information Sheet or contact us at techsupport@stemcell.com.

CARE OF MAGNET: After completing a cell separation, wipe the magnet clean with distilled water and then with alcohol to dry the magnet. To disinfect the magnet, wipe with 10% diluted household bleach and then wipe off with a water-dampened cloth to avoid corrosion of metal parts. Do not submerge the magnet in liquid.

PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED. FOR ADDITIONAL INFORMATION ON QUALITY AT STEMCELL, REFER TO [WWW.STEMCELL.COM/COMPLIANCE](http://www.stemcell.com/compliance).

Copyright © 2021 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and EasySep are trademarks of STEMCELL Technologies Canada Inc. Corning is a registered trademark of Corning Incorporated. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.