

AggreWell™ Rinsing Solution



Rinsing solution for AggreWell™ plates

Catalog # 07010

100 mL

Scientists Helping Scientists™ | WWW.STEMCELL.COM

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

Product Description

AggreWell™ Rinsing Solution is a surfactant solution that is required when using AggreWell™ plates to form embryoid bodies (EBs) or spheroids. Using AggreWell™ Rinsing Solution lowers the surface tension of AggreWell™ plates, making it easier to remove bubbles from the microwells. It also prevents cell adhesion to the plate and promotes efficient EB and spheroid formation. The lowered surface tension makes it easier to harvest EBs or spheroids from the microwells.

Properties

Storage: Store at 15 - 25°C.

Shelf Life: Stable until expiry date (EXP) on label.

Handling / Directions For Use

NOTE: AggreWell™ Rinsing Solution is required when using AggreWell™ plates.

1. Rinse AggreWell™ plates by adding AggreWell™ Rinsing Solution to each well to be used, as follows:
 - For 24-well plates: 0.5 mL/well
 - For 6-well plates: 2 mL/well
2. Centrifuge at 2000 x g for 5 minutes to displace bubbles.
3. Remove AggreWell™ Rinsing Solution from wells.
4. Wash each well to be used with 2 mL of DMEM/F-12 or phosphate-buffered saline (PBS) prior to adding cells.

For complete instructions for using AggreWell™ Rinsing Solution, refer to the Product Information Sheet for AggreWell™800 (Document #DX21397), available on our website at www.stemcell.com or contact us to request a copy.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2016 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and AggreWell are trademarks of STEMCELL Technologies Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.