

# Horse Serum for Human Myeloid Long-Term Cultures

For preparation of human long-term culture medium

Catalog # 06750

500 mL



Scientists Helping Scientists™ | [WWW.STEMCELL.COM](http://WWW.STEMCELL.COM)

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

[INFO@STEMCELL.COM](mailto:INFO@STEMCELL.COM) • [TECHSUPPORT@STEMCELL.COM](mailto:TECHSUPPORT@STEMCELL.COM)

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

## Product Description

This horse serum has been tested with all the reagents used in the preparation of MyeloCult™ H5100 (Catalog #05100). Components combined with this horse serum should be pre-tested in your culture system and selected to provide optimal support of human hematopoietic progenitors in long-term cultures.

This product is recommended for use in myeloid long-term culture media for human cells.

## Properties

**Storage:** Store at -20°C.

**Shelf Life:** Not applicable

## Handling / Directions For Use

1. Thaw at room temperature (15 - 25°C) or overnight at 2 - 8°C.
2. Swirl the bottle to mix.
3. Store at 2 - 8°C for up to 2 months. Alternatively, aliquot and store at -20°C. After thawing aliquots, do not re-freeze.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2016 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and MyeloCult are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.