

Collagenase/Hyaluronidase



Scientists Helping Scientists™ | WWW.STEMCELL.COM

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

10X Collagenase/hyaluronidase in DMEM

Catalog # 07912

10 mL

Product Description

10X Collagenase/Hyaluronidase in Dulbecco's Modified Eagle's Medium (DMEM) for the enzymatic dissociation of human and mouse mammary tissue.

Properties

- Storage:** Store at -20°C.
- Shelf Life:** Stable until expiry date (EXP) on label.
- Contains:**
- 3000 U/mL Collagenase
 - 1000 U/mL Hyaluronidase
 - DMEM (1000 mg D-glucose/L)

Materials Required But Not Included

- Ammonium Chloride Solution (Catalog #07800)
- EpiCult™-B Medium Kit (Human) (Catalog #05601)
- EpiCult™-B Mouse Medium Kit (Catalog #05610)
- HBSS with 10 mM HEPES, Without Phenol Red (Catalog #37150)
- Tissue Dissociation Flask (Catalog #27300)

Handling / Directions For Use

Thaw Collagenase/Hyaluronidase at room temperature (15 - 25°C) or overnight at 2 - 8°C.

NOTE: Collagenase/Hyaluronidase is provided as a 10X stock solution, and should be diluted 1 in 10 as outlined in the protocol below. Once thawed, use immediately or aliquot and store at -20°C until the expiry date as indicated on the label. After thawing aliquots, use immediately. Do not re-freeze.

NOTE: Avoid the use of glass pipettes and tubes when handling mammary epithelial cells, as these cells will stick to glass.

DISSOCIATION OF HUMAN MAMMARY TISSUE

1. Transport human mammary tissue from the operating room on ice in sterile specimen cups in complete EpiCult™-B Medium (Human) supplemented with 5% fetal bovine serum (FBS).
2. Transfer tissue to sterile glass petri dishes, mince with scalpels and then transfer to tissue dissociation flasks.
NOTE: Glass petri dishes can be used for this initial dissociation, as the concentration of epithelial cells is very low.
3. Dilute 1 part Collagenase/Hyaluronidase (10X stock) with 9 parts complete EpiCult™-B Medium (Human) and add to the minced tissue in the dissociation flasks. Ensure that the tissue is well-suspended in the enzyme mixture and the final volume is level with the widest portion of the flask. Cover the opening of the flask with sterile aluminum foil.
4. Gently dissociate the minced tissue on a rotary shaker at 37°C until all large tissue fragments are digested. Typical digestion time is 16 hours (overnight) for normal human mammary tissue. Longer digestion times may be required for tough fibrous tissue, shorter digestion times for softer tissue.
NOTE: The flasks should be sealed if the rotary shaker is not in a 5% CO₂ incubator.
5. Transfer the dissociated tissue to 50 mL centrifuge tubes, and centrifuge at 80 x g for 30 seconds.
6. Discard the overlying liquefied fat layer. The pellet ("A" pellet) is highly enriched for epithelial organoids.
7. Transfer the supernatant to a new 50 mL centrifuge tube and centrifuge at 200 x g for 3 minutes. The pellet ("B" pellet) from this second centrifugation contains variable numbers of epithelial cells, stromal cells, and red blood cells.

8. The supernatant from the second centrifugation is enriched for human mammary fibroblasts. To collect, transfer the supernatant to a new 50 mL centrifuge tube and centrifuge at 350 x g for 5 minutes.
9. The different cell fractions can now be cryopreserved. It is recommended that cells are cryopreserved in complete EpiCult™-B Medium (Human) supplemented with 50% FBS and 6% dimethyl sulfoxide (DMSO).

To generate a single-cell suspension from the “A” or “B” pellet, refer to the Technical Bulletin: A Guide to Solid Tissue Dissociation (Document #29107), available on our website at www.stemcell.com, or contact us to request a copy.

DISSOCIATION OF MOUSE MAMMARY TISSUE

1. Dilute 1 part Collagenase/Hyaluronidase (10X stock) with 9 parts complete EpiCult™-B Medium (Mouse) supplemented with 5% FBS in a 14 mL or 50 mL tube.
NOTE: Approximately 2 - 5 mL of the EpiCult™-B Medium/Collagenase/Hyaluronidase/FBS solution will be required for every 2 mammary glands to be dissociated.
2. Resect mammary glands and transfer to a sterile glass petri dish. Mince with scalpels in a cross-wise pattern until glands are rendered to a paste. Transfer the mammary tissue to the tube containing EpiCult™-B Medium/Collagenase/Hyaluronidase/FBS and incubate at 37°C for 6 - 8 hours with occasional pipetting and vortexing.
3. Centrifuge the cells at 350 x g for 5 minutes and discard the supernatant.
4. Resuspend the cell pellet in a mixture of 1 part cold Hanks' Balanced Salt Solution (HBSS) with 10 mM HEPES, Without Phenol Red supplemented with 2% FBS and 4 parts Ammonium Chloride Solution and centrifuge at 350 x g for 5 minutes. The resultant pellet contains epithelial cell organoids as well as stromal cells and lymphocytes.

To generate a single-cell suspension of mammary epithelial cells, refer to the Technical Bulletin: A Guide to Solid Tissue Dissociation (Document #29107), available on our website at www.stemcell.com, or contact us to request a copy.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2016 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and EpiCult are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.